


Changing the bulb in the Dolan Jenner halogen light source


1. **** Lamp Housing can still be very hot! Before changing the bulb make sure the lamp has been turned off for at least 10 minutes!!** Loosen the 2 screws on the front of the lamp housing all of the way. Indicated here with the red arrows.

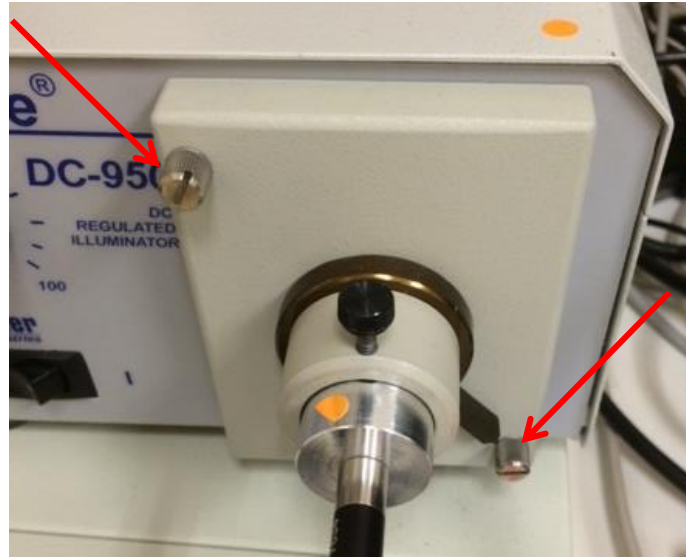


2. Pull the lamp housing out to expose the socket that connects to the bulb.



3. Slide the bulb out of its holder between the metal plate and 2 retaining clips.

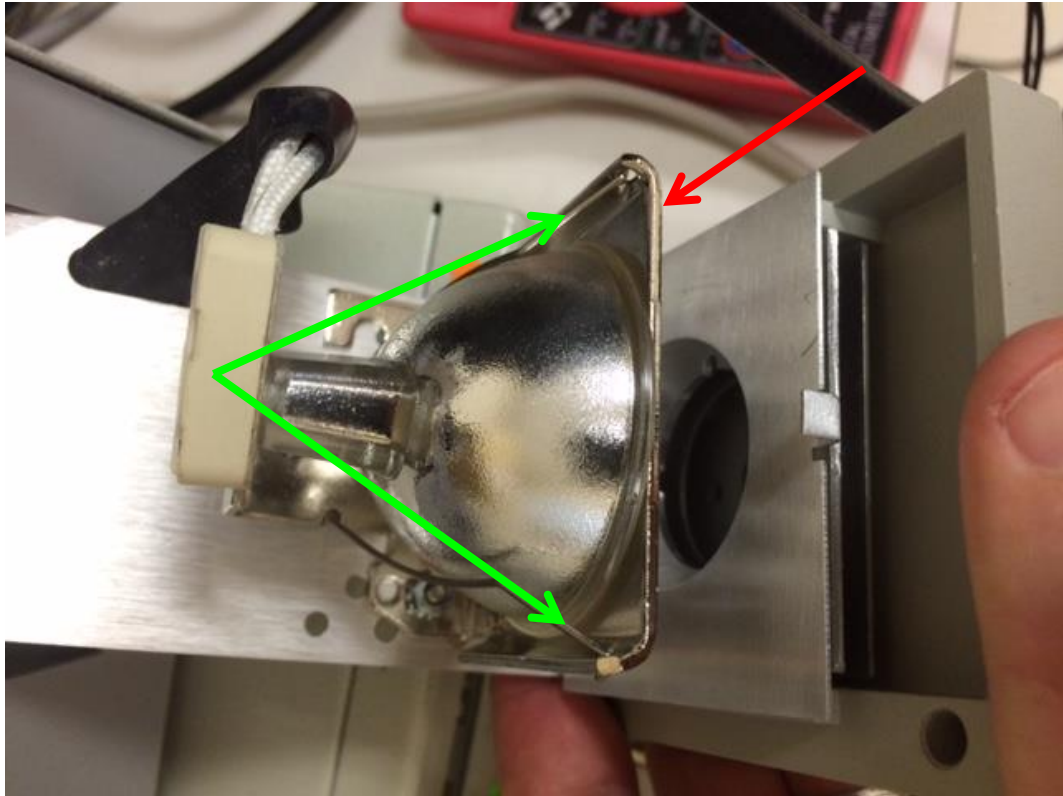


4. Gently pull the bulb from the socket. You may have to pull and rock the socket back and forth to set free.

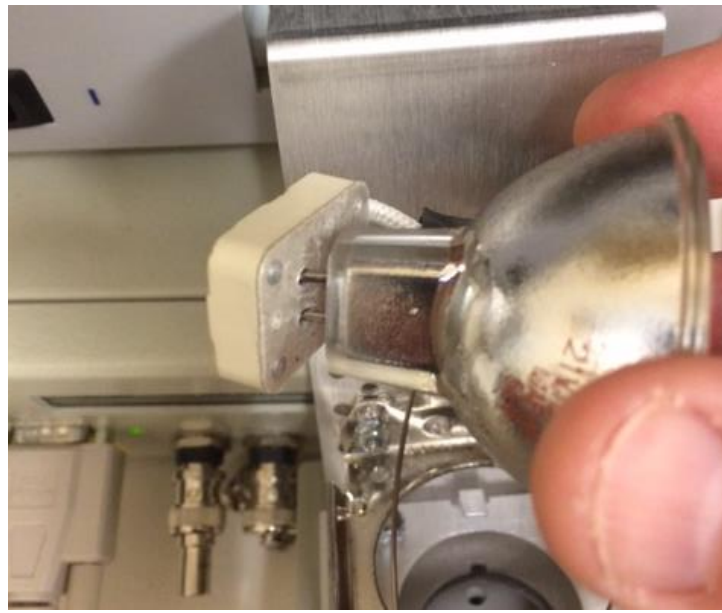

5. Place the new bulb in the socket and make sure the 2 prongs are all the way in the ceramic plug.



6. Place the bulb back in its housing sliding it between the metal plate and the 2 retaining clips.

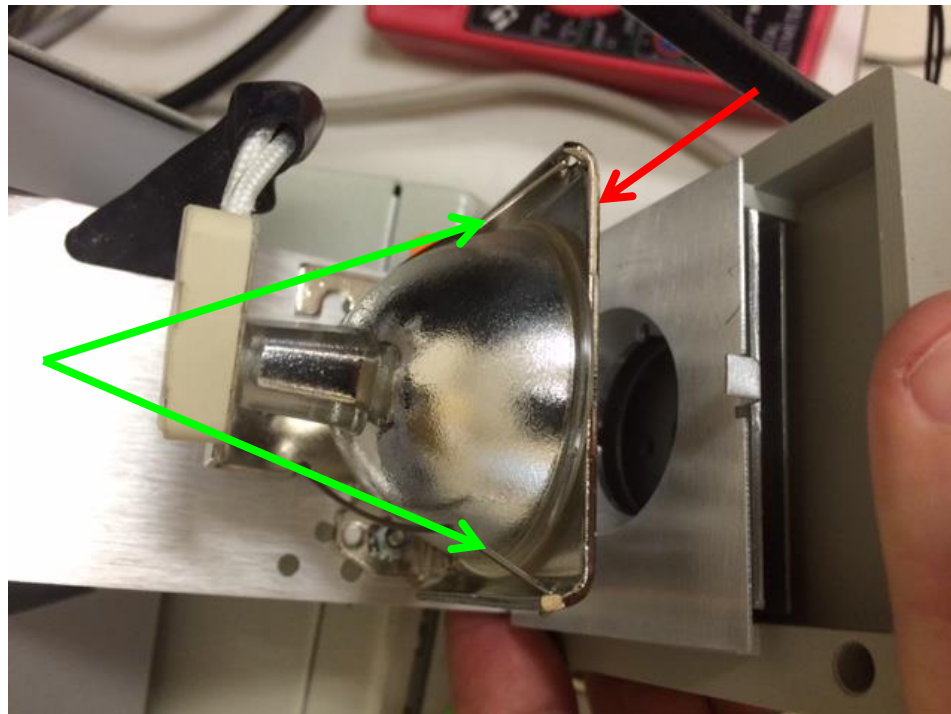

7. Push lamp housing back up against main body and tighten both screws.

